GridION

Quick start guide

Pre-install

This Quick start guide sets up your GridION™ for use.

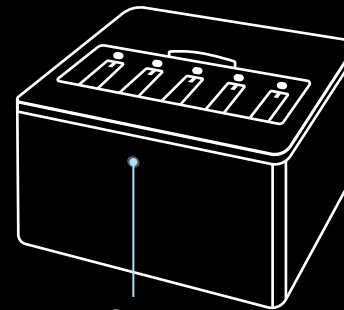
See the following:

GridION user manual
community.nanoporetech.com/to/gridion

Safety and regulatory information
community.nanoporetech.com/to/safety

For troubleshooting, see the user manual.

GridION

Power button

What's in the box:

GridION

Configuration Test Cells (x5)

Power cables*

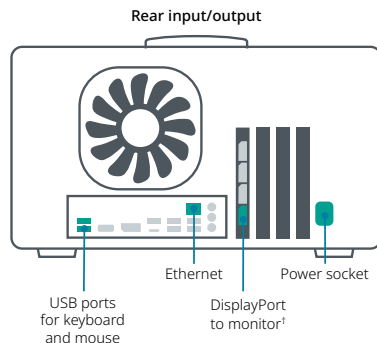
DisplayPort to HDMI adapter

*The GridION ships with 5 x power cables (1 US, 1 UK, 1 EU, 1 CN, 1 AUS) for international use.

1

Set up your device

1. **Unpack** your GridION device*.
2. Attach the **cables** and **peripherals** (as shown opposite).
3. Connect your power cable to your **power socket** and **power supply** and switch on the mains, as needed.
4. Press the **power button** (as shown on page 3).

*Ensure 30 cm clearance to front, rear and sides; keep away from direct heat and sunlight. See user manual for detailed advice.

¹If using HDMI only monitor, use the included DisplayPort to HDMI adapter.

! Only use the ports and connections coloured in green for installation.

2

Log in to MinKNOW

1. **Log in** to your GridION
Username: **grid**
Password: **grid**
2. Update your GridION **password**.
Your new password must:
 - Have a minimum 8 characters
 - Contain a capital letter and a lower-case letter
 - Contain a number or special character
 - Not be a palindrome (e.g. "radar")
 - Differ from a previous password by more than 2 characters.
3. Click  on the desktop to load MinKNOW™.
4. Log in to MinKNOW using your Oxford Nanopore account details.
5. Familiarise yourself with the tutorials (top right) or click **Skip all tutorials**.

3

Update software

1. Select **Device settings** in the side bar.
2. Click **Software** in the side bar.
3. Click **Install update** in the pop-up.
4. Shut down your GridION using the **power off** instructions (Power off step 4).

4

Power off

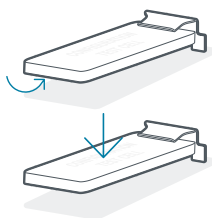
1. Select **Device settings** in the side bar.
2. Click **Shut down** in the main window.
3. Click **Shut down**  in the pop-up box.

Wait 10 seconds before turning on your GridION by pressing the power button. Repeat step 2 (Log in to MinKNOW) and then proceed to step 5 (Perform a Hardware check).

5

Perform a Hardware check

1. Slide the GridION lid back, insert the CTCs under each clip in positions X1-X5 (as shown) and slide the lid shut.

In MinKNOW:

2. Click **Start** (side bar).
3. Click **Hardware check** (under Utilities).
4. Click **Select all positions**.
5. Click **Start** (bottom right).
6. Confirm **Selected position(s) passed all checks** is displayed*.
7. Then, go to **Sequencing overview**.
8. Remove the CTC by reversing the movement to insert the CTCs.

Your GridION is now ready to start sequencing!

*If any hardware check fails: 1. Check the CTCs are fully inserted. 2. Repeat the Hardware check. See **Support** in **Additional information** if the problems persist.

6

The Nanopore Community

community.nanoporetech.com

Discover the Nanopore Community to find protocols, forums, support and downloads for your experiments.

TIP: Learn how to analyse your nanopore data at: nanoporetech.com/analyse

Additional information

- ✓ **Warranty**
A license and warranty can be purchased for your device here:
store.nanoporetech.com/device-warranty.html
Flow cell warranty, visit: community.nanoporetech.com/to/warranty
- 🔄 **Recycle used flow cells**
Oxford Nanopore is committed to environmental sustainability. You can help by sending your flow cells for recycling. Find out how: community.nanoporetech.com/support/returns
- 🛒 **Place your next order**
Buy more consumables at the Oxford Nanopore Store:
store.nanoporetech.com
- 📖 **Documentation**
Documentation for your device is available on the Nanopore Community:
community.nanoporetech.com/docs
- 👍 **Support**
For all of your customer and technical support needs, visit:
community.nanoporetech.com/support

Technical specification

Model number	GRD-MK1	
Supply voltage (V)	100-240 AC ± 10% (50-60 Hz)	
Maximum rated current (A)	6.5	
Maximum rated power (W)	650	
Size (H x W x D) (mm)	220 x 365 x 370	
Weight (kg)	14.4	
Installation ports	1 x Power socket 1 x Ethernet port (1 Gbps) 1 x Display port (for monitor)	1 x USB-A 2.0 port (for mouse) 1 x USB-A 2.0 port (for keyboard)
Software installed	Ubuntu, GridION OS, MinKNOW	
Compute specification	7 TB SSD storage, 64 GB RAM, Intel CPU (8-cores or more), 1 x NVIDIA GPU	
Environmental conditions	Functional range of electronics is within environmental temperatures of +5°C to +40°C. Users should allow 30 cm clearance to the rear and sides of the device. Designed to sequence in environmental temperatures of +18°C to +25°C. Intended for indoor use. Can be used up to altitudes of 2,000 m. Use within 30%–75% relative non-condensing humidity limits. The device has a Pollution Degree 2. WARNING: Rear of instrument heats up during operation.	

Oxford Nanopore Technologies

phone +44 (0)845 034 7900

email support@nanoporetech.com

 oxford-nanopore-technologies

 @nanopore

 @nanoporetech.com

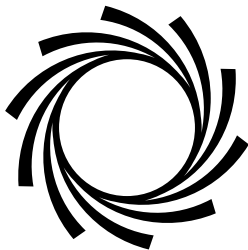
Gosling Building,
Edmund Halley Road,
Oxford Science Park,
Oxfordshire,
OX4 4DQ, UK

www.nanoporetech.com

Oxford Nanopore Technologies, the Wheel icon, GridION, and MinKNOW are registered trademarks or the subject of trademark applications of Oxford Nanopore Technologies plc in various countries. Information contained herein may be protected by patents or patents pending of Oxford Nanopore Technologies plc. All other brands and names contained are the property of their respective owners. © 2025 Oxford Nanopore Technologies plc. All rights reserved. Oxford Nanopore Technologies products are not intended for use for health assessment or to diagnose, treat, mitigate, cure, or prevent any disease or condition.

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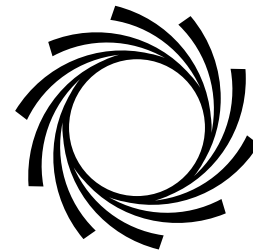
community.nanoporetech.com/to/gridion

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community.nanoporetech.com/to/safety

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